

GUÍA DE INICIO RÁPIDO

IMPRESIÓN Y ENVÍO DE INFORMES POR CORREO ELECTRÓNICO

Aprenda cómo:

- **1.** Para imprimir y descargar varios informes
- 2. Para enviar informes por correo electrónico



C LA	astrol BCHECK Fertung Labeheck Intelligence			Need Help	p 🕐 Search	😩 Acme Demo 👻
HOM	IE ACTIONS SAMPLES LABELS E	QUIPMENT ANALYTICS REPORTS DIS	TRIBUTION ADMIN PREFERENCE	S		
+	Add Filter					T
DATE	RANGE Last Year				🔀 EXPORT XLS 🚺 PRINT R	REPORT(S)
ļ	LAB NUMBER	SEVERITY ÷ C		CONDITION	UNIT ID	COMPONENT
	202309080644	Α	Acme Demo	Acceptable	221	ENGINE
	202309081234	Α	ACME Demo	Acceptable	293	ENGINE
	202309080603	C	ACME Demo	Wear	3093	HYDRAULIC
	202309081231	Α	ACME Demo	Acceptable	3021	ENGINE
	202309081233	Α	ACME Demo	Acceptable	297	ENGINE
	202309080611	D	ACME Demo	Viscosity/Wear	211	HYDRAULIC
	202309080641	C	ACME	Wear	211	ENGINE
	202308300905	B	ACME Dem		informes deada la	

Para imprimir uno o varios informes desde la pestaña MUESTRAS, seleccione las casillas de verificación situadas junto a las muestras y haga clic en el botón *IMPRIMIR INFORME(S)*.



															12	
E Ca	astrol					Bure	au Veril	tas Oil C	ondition N	lonitoring			San	ple A	nalysis	Repor
Account Information				1	2715 R	loyal Driv	/e, Staffor	d, TX					St	atus		
				lah	bocks	800-240	3-1118	itos com					Comple C			
				lau	Checksu	thhour	Dureauver	las.com					Sample C	ondition		
						Sar	nple Informa	tion			0	ther Sample	Information	1		
					-	Lab No.:	2019072906	12				PO No.:				
	Company Name Worksite	E ACME Del	TY				Sample 1	fracking #:	E201907180	592			Work Ore Referen	ler No.:	1028	
	Address	5: 123,	17				Rece	ived Date:	07/29/2019				Filt	er Age: 0	4020	
		Edmonton	, Alberta, Té	6B 3M9			Comple	eted Date:	08/01/2019				Make	Up Oil 0		
		nit Informat	ion					Com	anant Infor	nation			1	mount:	restion	
	Unit ID	nit mornat		ACTINAN		Com	nament D	com		ACTEM			Chuid Manuf		VDON	
	Unit Manufacture	r: Caterpillar	ERFUDDE	AST/MAI	IN	Comp	ponent Mar	nufacturer:	Caterpillar	(SIEM)			Fluid Manui	Fluid HEA	VRON VY DUTY PF	ANTIFR
	Unit Mode	1: 3516				2022010-2	Compon	ent Model:	3516				Brand/F	Product:		
	Unit Seria	I: UnivC					Compon	ent Serial:	UnivC	1223217			Fluid	Grade:		
	Unit WURShe	: Houston ,	IX				Compo	nent Type.	COOLING 3	SIEM						
Evaluate	d By: ENELSON	of S	AMING, whi SCA. Resam	ich can ca nple at nor MILLION	ause loss o rmal interv	f heat transfer al.	if this leve	el increase	s, Note a cha	nge in the co	olant color. 1	The coolant a	ppears stable.	Suggest add	ding full maint	GHT enance dose
Evaluate SPECTRO	d By: ENELSON	of S	AMING, whi SCA. Resam ARTS PER I	ich can ca nple at nor MILLION	ause loss o rmal interv	f heat transfer al.	if this leve	el increase	s, Note a cha	nge in the co	olant color. 1	The coolant a	ion protection ppears stable.	Suggest add	Jw. Note SLI ling full maint	GHT enance dose
Evaluate SPECTR LAB NO.	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN		AMING, whi CCA. Resam	ich can ca nple at nor MILLION	use loss o rmal interv	f heat transfer al.	if this levi	uce is rule	Ele	ments	unit color. 1	The coolant a	ppears stable.	Suggest add	un note SLI	GHT enance dose
LAB NO.	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019		AMING, whi CA. Resam	ich can ca nple at nor MILLION	rmal interv	of heat transfer al.	راب المراجع (f this level) رواب المراجع (f this level) رواب المراجع (f this level)	lincrease	S, Note a cha	ments	unit color. T	The coolant a	son protection ppears stable.	Suggest add	VV. Note SLI	GHT enance dose
LAB NO. 0612 1081	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019		AMING, whi SCA. Resam	ICh can ca nple at nor MILLION	nuse loss o rmal interv	theat transfer al.	ی پرچ <1	lincrease	S, Note a cha	ments	spinorous priorous pr	The coolant a	stable.	Suggest add	VV. Note SLI ling full maint	CHT enance dose
LAB NO. 0612 1081 0741	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019		AMING, whi SCA. Resam	ich can ca nple at nor MILLION	ause loss o rmal interv	theat transfer al.	ی پ ج ا	60000000000000000000000000000000000000	S, Note a cha	ments 1619 1659 1747	spinous provides prov	207 187 226	stable.	Suggest add	VV. Note SLI ling full maint	CHT enance dose
LAB NO. 0612 1081 0741 0820	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019	+0.0 of S LLYSIS IN PA & & & <1 <1 <1 <1 <1	AMING, whi SCA. Resam	ch can ca nple at nor MILLION S C C C C C C C C C C C C C C C C C C	ause loss o rmal interv	theat transfer al.	ی پ چ دا	6 6 7 7 16 15 17 17	Ele 6 334 335 359 360	ments 1619 1659 1747 1752	58 58 58 55 59	207 187 226 228	son protection ppears stable.	Suggest add √ √ <1 <1 <1 1 <1	VV. Note SLI ling full maint	CHT enance dose
LAB NO. 0612 1081 0741 0820 0873	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019	+0.0 of S LLYSIS IN PA <1 <1 <1 <1 <1 <1 <1	AMING, whi SCA. Resam	ch can ca nple at nor MILLION S C C C C C C C C C C C C C C C C C C	ause loss o rmal interv 1 1 1 1 1 1 1	theat transfer al.	ی پ چ دا	6 6 7 16 15 17 17 16	Elc	ments 1619 1659 1747 1752 1536	unit color. 1	207 187 226 228 204	65 68 67 66 61	Suggest add √ <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	<pre>VV. Note SLI ling full maint vio vio vio vio vio vio vio vio vio vi</pre>	CHT enance dose
LAB NO. 0612 1081 0741 0820 0873 0542	SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019	+0.0 of S LYSIS IN PA <1 <1 <1 <1 <1 <1 <1 <1	AMING, whi GCA. Resam ARTS PER 1 4000 4000 4000 4000 4000 4000 4000 40	ch can ca nple at nor MILLION	ause loss o rmal interv	theat transfer al.	ی پ درا	6 6 7 7 16 15 17 17 16 16	Elc	ments 1619 1659 1747 1752 1536 1676	58 58 58 65 59 42 52	207 187 226 228 204 218	100 protection ppears stable. 00 00 00 00 00 00 00 00 00 00 00 00 00	Suggest add √ √1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	Constraints of the second seco	
LAB NO. 0612 1081 0741 0820 0873 0542	SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 INFORMATION	FO. of S LYSIS IN PA S S N PA S S N PA S S N PA S S S N PA S S S S S S S S S S S S S S S S S S	AMING, whi GCA. Resam ARTS PER 1 4000000000000000000000000000000000000	ich can ca nple at nor MILLION	ause loss o rrmal interv	theat transfer al.	ی پ دا FLUID	6 6 7 7 16 16 15 17 17 16 16 16 0 PROPER	Ele	ments	(1) (1) (1) (1) (1) (1) (1) (1) (1) (1)	207 187 226 228 204 218	10n protection ppears stable. 00 00 00 00 00 00 00 00 00 00 00 00 00	vers are co Suggest add √ √ 1 <1 1 <1 <1 <1 <1 <1 <1 <1	Constraints of the second seco	GHT enance dose <1
LAB NO. 0612 1081 0741 0820 0873 0542 AMPLE Lab	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 INFORMATION Sample Sample	+0. of S LYSIS IN PA <1 <1 <1 <1 <1 <1 <1 <1 Unit	AMING, whi GCA. Resam ARTS PER 1 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4	In the second se	Filter	theat transfer al.	ی ج دا دا دا	6 16 15 17 16 16 15 17 16 16 16 17 17 16 16 16 17 17 16 16 16 17 17 16 16 16 17 17 17 16 16 16 17 17 17 16 16 17 17 16 16 17 17 17 16 16 17 17 17 16 16 17 17 17 16 16 17 17 17 16 17 17 17 16 17 17 17 17 17 17 17 17 17 17 17 17 17	Elc	ments 1619 1659 1747 1752 1536 1676	58 58 58 65 59 42 52	207 187 226 228 204 218	100 protection ppears stable. 000 000 000 000 000 000 000 000 000 0	Suggest add √ √ √ 1 <1 1 <1 <1 <1 <1 <1 <1 <1	VV. Note SLI ling full maint	Image: Constraint of the second se
LAB NO. 0612 1081 0741 0820 0873 0542 AMPLE Lab No. 0612	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 02/27/2019 INFORMATION Sample Drawn 07/10/2019	+0. of S LYSIS IN PA <1 <1 <1 <1 <1 <1 <1 <1 Unit Time 34829	AMING, whi GCA. Resam ARTS PER 1 4000 4000 4000 4000 4000 4000 4000 40	In the second se	Filter Chgd. No	theat transfer al.	ی ج ا	6 6 7 7 7 7 7 7 6 16 15 17 17 16 16 16 9 PROPER	Elc	ments	58 58 58 59 42 52	207 227 228 204 218	00 protection ppears stable. 00 00 00 00 00 00 00 00 00 00 00 00 00	Suggest add Suggest add √ √ <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	VV. Note SLI ling full maint	Image: Constraint of the second se
LAB NO. 0612 1081 0741 0820 0873 0542 AMPLE Lab No. 0612 1081	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 02/27/2019 02/27/2019 INFORMATION Sample Drawn 07/10/2019 06/28/2019	FO. of S LYSIS IN PA <1 <1 <1 <1 <1 <1 <1 Unit Time 34829 34541	AMING, whi GCA. Resam ARTS PER 1 ARTS PER 1 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4 4	In the second se	Filter Chgd. No	theat transfer al.	ی ج ا ا ا ا ا ا ا ا ا ا ا ا ا ا ا ا ا ا	6 6 7 7 7 7 7 6 6 7 7 7 7 7 7 7 6 7	Elc	ments 1619 1659 1747 1752 1536 1676 1676 1676	sprint color. 1 olant color. 1 58 58 58 59 42 52 42 52	207 187 226 228 204 218	65 68 67 66 61 64	Suggest add Suggest add <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	VV. Note SLI ling full maint vite	enance dose
LAB NO. 0612 1081 0741 0820 0873 0542 AMPLE Lab No. 0612 1081 0612 1081	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 04/11/2019 04/11/2019 04/12/2019 02/27/2019 INFORMATION Sample Drawn 07/10/2019 06/28/2019 06/28/2019 06/28/2019	CLYSIS IN PA	AMING, whi GCA. Resam ARTS PER 1 ARTS PER 1	In the second se	Filter Chgd. No No	<pre> theat transfer al. theat transfer al.</pre>	ج د دا الالالالالالالالالالالالالالالالال	00021514 el increase 16 15 17 17 16 16 16 16 16 0 PROPER	Elc	ments Market Solution ments Market Solution ments	olant color. T	207 187 226 204 218	65 68 67 66 61 64	<pre> Suggest add Suggest add cl cl</pre>	Construction of the second sec	enance dose
Valuate PECTR NO. 0612 1081 0741 0820 0873 0542 AMPLE Lab No. 0612 1081 0741 0820	d By: ENELSON OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 05/17/2019 04/11/2019 04/11/2019 04/27/2019 03/25/2019 02/27/2019 INFORMATION Sample Drawn 07/10/2019 06/28/2019 06/28/2019 05/17/2019	Clysis in PA Clysis in PA Cl	AMING, whi GCA. Resam ARTS PER I	In the second se	Filter Chgd. No No No	theat transfer al. theat transfe	ی ج دا ا	6 6 7 7 7 7 7 7 6 6 7 7 7 7 7 7 7 7 7 7	Solution of the second	ments 1619 1619 1659 1747 1752 1536 1676 1676	58 58 58 65 59 42 52	207 187 226 228 204 218	65 66 67 66 61 64	<pre> Suggest add Suggest add cl cl</pre>	Construction of the second sec	<pre>clipselogged clipselogged</pre>
LAB NO. 0612 1081 0741 0820 0542 AMPLE Lab No. 0612 1081 0741 0810 0741 0820 0873	SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 03/25/2019 03/25/2019 03/25/2019 03/25/2019 03/25/2019 03/25/2019 07/10/2019 06/28/2019 05/17/2019 06/28/2019 05/17/2019 06/28/2019 05/17/2019 04/11/2019<	CLYSIS IN PA	AMING, whi CCA. Resam ARTS PER 1 ARTS PER 1	In the second se	Filter Chgd. No No No No No	theat transfer al. theat transfer transfer al. theat transfer al. theat transfer transfer transfer transfer	یک دا ا	e ab	s, Note a cha Ele 334 335 359 360 313 341 TIES TITÁ U mprin	ments 1619 1619 1659 1747 1752 1536 1676 1676 1676	58 58 58 59 42 52 IEVA	207 187 226 228 204 218 0esta 0ne e	65 68 67 66 61 64 Ma coi	<pre> Suggest add Suggest add i i i i i i c f el P o de </pre>	vv. Note SLI ling full maint state of the state of the s	<pre>GHT enance dose</pre>

letra P.

KEY: UoM - Unit of Measure Y - Yes N - No C - Changed S - Sampled >

This analysis is intended as an aid in predicting mechanical wear. Test results, mai and apply only to this sample as provided. No guarantee, expressed or implied, is equipment and all of its components is the responsibility of the equipment owner. Testing performed by Bureau Veritas, an ISO/IEC 17025/2017 accredited laboratory by A accreditation. Click here for Tests and Methodologies

Para descargar, seleccione el icono de descarga.

N.



	Castrol ABCHECK Fortung Labebeck Intelligence				Need Help ⑦ Search	😩 Acme Demo 👻
HO	ME ACTIONS SAMPLES LABELS	EQUIPMENT ANALYTICS REPORT	s distribution admin prefe	RENCES		
+	Add Filter				1	T
DATI	E RANGE Last Year SHOW		/		EXPORT XLS	PRINT REPORT(S)
l		SEVERITY - Select -	CUSTOMER	CONDITION		
	202309080644	A	Acme Demo	Acceptable	221	ENGINE
	202309081234	A	ACME Demo	Acceptable	293	ENGINE
0	202309080603	c	ACME Demo	Wear	3093	HYDRAULIC
	202309081231	Α	ACME Demo	Acceptable	3021	ENGINE
	202309081233	Α	ACME Demo	Acceptable	297	ENGINE
\Box	202309080611	D	ACME Demo	Viscosity/Wear	211	HYDRAULIC
0	202309080641	С	ACME	Wear	211	ENGINE
	202308300905	B	_			ENGINE

Para *Exportar* una o más muestras a un archivo Excel desde la pestaña MUESTRAS, seleccione las casillas de verificación situadas junto a las muestras y haga clic en el botón *EXPORTAR XLS*.



e LA	astrol BCHECK Fortung Laboheck Intelligence				Need Help ⑦	Search	3 Downloading Report X
ном	IE ACTIONS SAMPLES LABELS E	QUIPMENT ANALYTICS REPORTS I	DISTRIBUTION ADMIN PREFEREN	ICES			
+	Add Filter						T
DATE	RANGE Last Year • SHOW				X	EXPORT XLS	PRINT REPORT(S)
0	LAB NUMBER	SEVERITY ÷		CONDITION - Select -	÷ UNIT	ID	COMPONENT
	202309080644	Α	Acme Demo	Acceptable	221		ENGINE
	202309081234	Α	ACME Demo	Acceptable	293		ENGINE
	202309080603	C	ACME Demo	Wear	3093		HYDRAULIC
	202309081231	Α	ACME Demo	Acceptable	3021		ENGINE
	202309081233	Α	ACME Demo	Acceptable	297		ENGINE
	202309080611	D	ACME Demo	Viscosity/Wear	211		HYDRAULIC
	202309080641	C	ACME	Wear	211		ENGINE
	202308300905	B	ACME Demo	Low Viscosity	202		ENGINE

El archivo Excel se creará y se descargará automáticamente; el enlace al archivo aparecerá en la parte inferior o superior de su página, dependiendo de su navegador.





RECOMMENDATIONS

Maintenance for Lab No.: 201907290612 Evaluated By: ENELSON ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED! Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of heat transfer if this level increases. Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose of SCA. Resample at normal interval.

> Para imprimir o descargar un informe desde la pantalla Detalles, pulse el botón **DESCARGAR PDF**.





RECOMMENDATIONS

Maintenance for Lab No.: 201907290612 Evaluated By: ENELSON ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED! Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of heat transfer if this level increases. Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose of SCA. Resample at normal interval.

> Si esta opción está seleccionada, se incluirán en el PDF los Gráficos, Anexos, Comentarios y Acciones. Su selección se conservará para futuras impresiones.



Control Contro Control Control Control Control Control Control Control Control Co	Construction Bureau Veritas Cil Condition Monitoring 12715 Royal Drive, Staffrid, TX 800-248-7778 Sample Analysis Report Status Construction Construction Construction Sample Analysis Report Status Construction Sample Analysis Report Status Construction Lab Customer (Dr. 30382 Construction Construction Construction Construction Market (Dr. 30382) Construction Construction Construction Construction Market (Dr. 30382) Construction Construction Construction Construction Construction Market (Dr. 30382) Construction Component Information Full Information Full Information Full Information Market (Dr. 1007CL/HER FLODE ASTIMALI) Unit Mandeture: Component Information Full Information Full Information Unit Mandeture: Component Information Component Information Full Information Full Information Market (Dr. 1007CL/HER FLODE ASTIMALI) Unit Mandeture: Component Information Full Information Full Information Market (Dr. 1007CL/HER FLODE ASTIMALI) Unit Mandeture: Component Information Full Information Full Information Market (Dr. 1007CL/HER FLODE ASTIMALI) Unit Mandeture: Component Serial Unit Construction Status Component Serial Unit Construction Status Component Network Status																	
Matchecksupport@bureauveritas.com Sample formation Account Information Company Name: ACME Eering Company Name: ACME Eering Company Nam	12715 Royal Drive, Stafford, TX 12715 Royal Drive, Stafford, TX 12715 Royal Drive, Stafford, TX 1280-248-7778 1280-248-7778 1280-248-7778 1280-248-7778 1280-248-7778 1280-248-7778 1280-248-7778 1280-228-7778 1280-228-7778 1280-228-7778 1280-228-2778 1280-228-2789 1280-228-2789 1280-228-2789 1280-228-2789 1280-228-2789 1280-228-2789 1280-228 1280-228 1280-228 1280-228 1280-228 1280-228 1280-228 1280-228 1280-228 1280-228 1280-228 1280-228 1280-2280 1280-2280 1280-2280 1280-2280 1280-2280 1280-2280 <th colspan<="" td=""><td>CC</td><td>astrol</td><td></td><td></td><td></td><td></td><td>Bureau Ve</td><td>ritas Oil C</td><td>ondition N</td><td>lonitoring</td><td></td><td></td><td>San</td><td>nple A</td><td>nalysis</td><td>Repor</td></th>	<td>CC</td> <td>astrol</td> <td></td> <td></td> <td></td> <td></td> <td>Bureau Ve</td> <td>ritas Oil C</td> <td>ondition N</td> <td>lonitoring</td> <td></td> <td></td> <td>San</td> <td>nple A</td> <td>nalysis</td> <td>Repor</td>	CC	astrol					Bureau Ve	ritas Oil C	ondition N	lonitoring			San	nple A	nalysis	Repor
Differentiation Sample Information Company home: Schedering Company home: Schedering Schedering <th col<="" td=""><td>Instruction Sample formation Account Information Sample Condition Account Information Sample Information Other Sample Information Description Sample Information Other Sample Information Policity Information Description Sample Information Other Sample Information Policity Information Description Sample Information Component Description: Status Information Fluid Manufacture: Clear Note Status Information Description Sample Information Component Description: Status Information Policity Information Description Status Information Component Description: Status Information Policity Information Policity Information Description Status Information Status Information Policity Information Policity Information Description Status Information Status Informat</td><td></td><td>RCHE</td><td>CK</td><td></td><td></td><td></td><td>12715</td><td>Royal Driv</td><td>e, Staffor</td><td>d, TX</td><td></td><td></td><td></td><td></td><td>SI</td><td>tatus C</td></th>	<td>Instruction Sample formation Account Information Sample Condition Account Information Sample Information Other Sample Information Description Sample Information Other Sample Information Policity Information Description Sample Information Other Sample Information Policity Information Description Sample Information Component Description: Status Information Fluid Manufacture: Clear Note Status Information Description Sample Information Component Description: Status Information Policity Information Description Status Information Component Description: Status Information Policity Information Policity Information Description Status Information Status Information Policity Information Policity Information Description Status Information Status Informat</td> <td></td> <td>RCHE</td> <td>CK</td> <td></td> <td></td> <td></td> <td>12715</td> <td>Royal Driv</td> <td>e, Staffor</td> <td>d, TX</td> <td></td> <td></td> <td></td> <td></td> <td>SI</td> <td>tatus C</td>	Instruction Sample formation Account Information Sample Condition Account Information Sample Information Other Sample Information Description Sample Information Other Sample Information Policity Information Description Sample Information Other Sample Information Policity Information Description Sample Information Component Description: Status Information Fluid Manufacture: Clear Note Status Information Description Sample Information Component Description: Status Information Policity Information Description Status Information Component Description: Status Information Policity Information Policity Information Description Status Information Status Information Policity Information Policity Information Description Status Information Status Informat		RCHE	CK				12715	Royal Driv	e, Staffor	d, TX					SI	tatus C
Account information Sample Information Other Sample Information Liab No: 201907200512 PO No: Worksite: Houston, TX Sample Tarking #: Ecologit20150592 Sample Tarking #: Ecologit20150592 Worksite: Houston, TX Sample Tarking #: Ecologit20150592 Sample Tarking #: Ecologit20150592 Worksite: Houston, TX Sample Tarking #: Ecologit20150592 Sample Tarking #: Ecologit20150592 Work Order Edmonton, Aberta, T6B 3M9 Component Information Fluid Monderson Unit Information Component Secretion: Component Information Fluid Monderson Unit Note: State Fluid Note: State Component Secretion: Component Mode: State Pluid Monderson Unit Note: State Inductore: Component Mode: State Fluid Monderson Unit Note: State Component Mode: State Pluid Monderson State Secretion: Component Mode: State Pluid Monderson	Account Information Sample Information Other Sample Information Lab No: 201907290512 Lab No: 201907290512 Provide	Featuring Labcheck Intelligence				labcheck	support@	oureauver	itas.com					Sample C	ondition			
Link Octoment Dis: 31912 Company Name: CALLE Damin Address: 123, Edmointon, Aberta, T6B 3M9 Lab No.: 301907280612 Sample Tracking #: E01907100592 Breceived Date: 0729/2019 Completed Date: 0729/2019 Component Model: 2516 Unit Model: 2516 Unit Serial UnivC Unit Model: 2516 Unit Serial UnivC Unit Model: 2516 Unit Serial UnivC Unit Model: 2516 File: Age: 0 Marker Up Oil 0 Amount: Marker Up Oil Unit Identification: Component Information Fluid Information Fluid Information Multi Manufacture: Categorian Unit Model: 2516 Unit Serial UnivC Unit Model: 2516 Component Model: 2516 Component Model: 2516 Fluid Information Internance for Lab No: 307220612 AMAU/SIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED/ Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of the internant interval. Fluid Manufacture: Categorian Fluid Information Internance for Lab No: 307220612 SCA. Resample at normal interval. Erements Erements Internance for Lab No: 307220612 GCA. Resample at normal interval. Erements Erements Internance for Lab No: 307220612 GCA. Resample at normal interval. Erements Erements Internance for Lab No: 307220612 GCA. Resample at normal interval. Erements Erements Internance for Lab No: 30722051 GCA. Resample at normal interval. Erements Erements	Lub Custom Company Name: 4/LINE Demoty Company Name: 4/LINE Demoty Address: 123, Edmotion, Aberta, TBB 3M9 Lub No: 2019/07/30632 Sample Tracking N = 2019/7130639 Received Date: 1/729/2019 Component Mode: 2019/20199 Received Date: 1/729/2019 Component Mode: S108 Unit Noted: S108 Unit Noted	Account Information					Sar	nle Informa	tion			0	ther Sample	Information				
Company Name: 40:MED terms Sample Tracking # E03907189592 Work Order No:: Reference No:: 758028 Work Name: Edmontion, Alberta, T6B 3M9 Sample Tracking # E03907189592 Sample Abs: 071292019 Sample Abs: 071292019 Unit Information Completed Date: 071292019 Sample Tracking # E03907189592 Sample Tracking # E03907189592 Sample Tracking # E03907189592 Unit Information Completed Date: 071292019 Sample Tracking # E03907189592 Sample Tracking # E03907189592 Unit Manufacture: Caterplitar Component Information Fluid Information Unit Manufacture: Caterplitar Component Mode: 3516 Component Mode: 3516 Unit Seria: Unit Seria: Component Mode: 3516 Component Mode: 3516 Fluid Alformation S0228021 Market Manufacture: Caterplitar Component Mode: 3516 Component Mode: 3516 Fluid Rand S02280221 Control Mode: SCA. Resample at normal interval. Component Mode: 3516 Component Mode: 3516 Component Mode: 3516 Component Mode: 3516 S02280221 Control Mode: Sca. Resample at normal interval. Component Mode: 3516 Component Mode: 3516 Component Mode: 3516 Component Mode: 3516 Component Mode: 3516 <td>Company Name: ACMEDiane Sample Tracking # E201907180992. Work Criter No: More No: Work Criter No: Received Date: 0779/02019 Received Date: 0779/02019 Work Criter No: Received Date: 0779/02019 Medica: Ethnomon, Alberta, T6B 3M9 Component Information Fluid Anulacture: Caterplilar Received Date: 0779/02019 Received Date: 0700/02019 Received Date: 0700/02019<!--</td--><td></td><td>Lab Customer ID#</td><td>319182</td><td>ation</td><td></td><td></td><td></td><td>Lah No.</td><td>2019072906</td><td>12</td><td></td><td></td><td></td><td>PO No .</td><td>monnation</td><td></td></td>	Company Name: ACMEDiane Sample Tracking # E201907180992. Work Criter No: More No: Work Criter No: Received Date: 0779/02019 Received Date: 0779/02019 Work Criter No: Received Date: 0779/02019 Medica: Ethnomon, Alberta, T6B 3M9 Component Information Fluid Anulacture: Caterplilar Received Date: 0779/02019 Received Date: 0700/02019 Received Date: 0700/02019 </td <td></td> <td>Lab Customer ID#</td> <td>319182</td> <td>ation</td> <td></td> <td></td> <td></td> <td>Lah No.</td> <td>2019072906</td> <td>12</td> <td></td> <td></td> <td></td> <td>PO No .</td> <td>monnation</td> <td></td>		Lab Customer ID#	319182	ation				Lah No.	2019072906	12				PO No .	monnation		
Worksite: Houston, TX Sampled Date: 07/10/2019 Reference No: 7594028 Adress: 123, Edmonton, Aberta, T6B 3M9 Component Information Filter Age: 0 Make Up OI 0 Amount Unit Information Component Information Fluid Information Unit ID: 1107/102109 Component Manufacture: Caterpline Unit Monté: 3316 Fluid Information Unit Manufacture: Caterpline Unit Monté: 3316 Component Manufacture: Caterpline Component Manufacture: Caterpline Unit Worksite: Hundon Fluid Manufacture: CHEVRON Unit Monté: 3316 Component Manufacture: Caterpline Component Manufacture: Caterpline Statepline Component Manufacture: Caterpline Statepline Component Manufacture: Caterpline Component	Worksite: Houston, TK Sampled Date: 07/10/2019 Reference No: 7594028 Bit Page 10 Sampled Date: 07/02/2019 Reference No: 7594028 Unit Hormation Component Information Full Harvan Unit Munitic: 1107/ELMER FUDD EASTMAIN Component Information Full Harvan Unit Munitic: 1107/ELMER FUDD EASTMAIN Component Manufacture: Caterplan Reference No: 7594028 Unit Munitic: 1107/ELMER FUDD EASTMAIN Component Manufacture: Caterplan Public Harvan Full Harvan Unit Model: 3516 Component Manufacture: Caterplan Component Manufacture: Caterplan Public Harvan Public Harvan Unit Model: 3516 Component Manufacture: Caterplan Component Manufacture: Caterplan Component Manufacture: Caterplan Public Harvan Unit Worksite: Houston, TX Marvan Component Manufacture: Caterplan Component Manufacture: Caterplan Public Harvan Unit Worksite: Nueston, TX Marvan Marvan Marvan Public Harvan Public Harvan 1907280612 Component Manufacture: Caterplan Marvan Public Harvan Public Harvan Public Harvan 1907280612 Component Manufacture: Caterplan Marvan Public Harvan Public Harvan		Company Name	ACME Der	mo			Sample	e Tracking #:	E201907180	592			Work Or	der No.:			
Address: 123, Edmontor, Alberta, T6B 3M9 Preceded Date: 07/29/0219 Print Age: 0 Make Up 01 0 Amount: Unit Information Component Information Filter Age: 0 Make Up 01 0 Amount: Print Age: 0 Make Up 01 0 Amount: Unit Moritature: Categoriest Date: 08/01/2019 Unit Moritatie: Component Information Filter Age: 0 Make Up 01 0 Amount: Unit Moritature: Categoriest Date: 08/01/2019 Unit Moritatie: Component Mandacture: Categoriest Component Markacture: Print Age: 0 Make Up 01 0 Amount: Print Age: 0 Make Up 01 0 Amount: Unit Moritatie: Statistic: Component Mandacture: Component Markacture: Component Markacture: Print Age: 0 Make Up 01 0 Amount: Print Age: 0 Make Up 01 0 Amount: Unit Moritatie: Statistic: Component Mandacture: Categoriest Categoriest Component Markacture: Statistic: Print Age: 0 Make Up 01 0 Amount: Print Age: 0 Age:	Address: Little Received Date: M22/93 Het Age: 0 Annout: Unit Information Component Information Fluid Manufacture: Amount: Unit Information: Component Information Fluid Manufacture: Fluid Manufacture: Unit Information: Component Description: COCULING SYSTEM Component Manufacture: Fluid Manufacture: Unit Model: 3516 Component Description: Component Thescription: Fluid Manufacture: Unit Model: 3516 Component Thescription: Component Thescription: Fluid Manufacture: Internance for Lab NC: DAMINOS, which can cause loss of heat transfer If this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance does of SCA Resample at normal interval. Internance for Lab NC: DAMINOS, which can cause loss of theat transfer If this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance does of SCA Resample at normal interval. EXTOCHEMICAL AMALYSIS IN PARTS PER MILLON Entertains AB SMARPE Age: 0 Age		Worksite	: Houston ,	ТХ			Sa	impled Date:	07/10/2019				Referen	nce No.: 759	4028		
Amount: Unit Mondature: Caterpillar Component Description: COOLING SYSTEM Unit Manufacture: Caterpillar Component Manufacture: Caterpillar Pluid Manufacture: CheVRON Unit Kode: 3516 Component Manufacture: Caterpillar Pluid Manufacture: CheVRON Fluid Manufacture: CheVRON Unit Kode: 3516 Component Serial: UnivC Component Monde Staff Fluid Manufacture: CheVRON Fluid Grade: Interance for Labor: ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED Mote some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT Fluid Manufacture: CheVRON Fluid Manufacture: CheVRON 100220512 Manufacture: Service Ser	Unit Manufacture: Component Information Fluid Information Unit D: 1107/ELMER FUDD EASTMAIN Component Manufacture: Camponent Manufacture: Cheven the state of the s		Address	Edmonton.	Alberta, T6E	3M9		Com	pleted Date:	07/29/2019 08/01/2019				Mak	ter Age: 0 e Up Oil 0			
Unit Information Component Information Fluid Information Unit Unit Dir Dir Dir LINGE LIND EASTMAIN Unit Manufacture: Caterpillar Unit Manufacture: Caterpillar Unit Modei: 3516 Unit Senia: Univol Unit Workite: Houston, TX Component Modei: 3516 Component Modei: 3516 Component Modei: 3516 Component Modei: 3516 Component Modei: 3516 Component Modei: 3516 Component Modei: 3516 Pluid Manufacture: Caterpillar Brand/Product: Fluid Grade: Unit Workite: Houston, TX ANLYSIS INDICATES COLANT MAINTENANCE/ SERVICE IS REQUIRED I Nete some phosphorous present. Corrosion protection levels are LOW. Note SLICHT FOAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dos of SCA. Resample at normal interval. CTCOCHEMICAL ANALYSIS IN PARTS PER MILLION Elements It 00/28/2019 c1 c1 c1 15 334 1619 58 207 65 c1 c	Unit Information Component Model: SDI Component Model: SDI Fluid Information Unit Wanufacture: Caterpillar Unit Model: SDI Unit Model: SDI Component Manufacture: Caterpillar Component Model: SDI Component Model: SDI Component Type: COOLING SYSTEM Fluid Manufacture: CHEVRON Fluid HEAVY DUTY PF ANTIFR Brand/Product: Fluid Grade: Internace for Lab No: B072290512 MALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED! Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT POAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose allose of SCA. Resample at normal interval. ECTROCHEMICAL ANALYSIS IN PARTS PER MILLION Elements Ass SAMPLE DRAWN add add add add add add allo 0628/2019 c1 c1 c1 15 335 1659 58 197 68 c1 c1 allo 0628/2019 c1 c1 c1 16 334 1619 58 207 65 c1 c1 c1 allo 0628/2019 c1 c1 c1 c1 16 334 1619 58 207 65 c1 c1 c1 allo 0628/2019 c1 c1 c1 c1 16 333 1356 42 204								•						Amount:			
Unit Bild 11077ELMEE PLOD EASTMAIN Component Description: COULING SYSTEM Fuld Manufacture: CHEVRON Unit Mode: 35.6 Component Mode: 35.16 Component Mode: 35.16 Bild Manufacture: CHEVRON Unit Mode: 35.6 Component Mode: 35.16 Component Mode: 35.16 Bild Manufacture: CHEVRON Unit Mode: 35.6 Component Mode: 35.16 Component Mode: 35.16 Bild Grade: Unit Worksite: Houston , TX ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED! Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FodMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dos 007290612 SAMPLE Seample at normal interval. Elements Elements State	Unit ID: ID:/ELMER PLODE EASTMAIN: Unit Mandiacture: Caterpillar Unit Mode: S516 Unit Serial: UnivC Unit Worksite: Houston, TX Component Description: COLUNG SYSTEM Fluid Mandiacture: CHEVRON ErandProduce: Unit Worksite: Houston, TX Internance for Lab No: Di07290512 AALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause bosh heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose alcaded By: ENELSON ECTROCHEMICAL ANALYSIS IN PARTS FER MILLION Efferentis ECTROCHEMICAL ANALYSIS IN PARTS FER MILLION Efferentis Ass SAMPLE & & & & & & & & & & & & & & & & & & &		Ui	nit Informati	ion				Comp	onent Inform	nation				Fluid Info	ormation		
Bit Mathematic Description Description Description Description Unit Model: 2516 Component Seria: Unit Drawn The analysis Drawn Dr	Component Material Component Material Component Material Component Material Component Material Component Serial: Dirac Brand/Product: Full Brand/Product: 10012 Unit Morkits: House Component Serial: Unit Component Serial: Dirac Di		Unit ID	: 1107/ELM	ER FUDD EA	ST/MAIN		Component N	Description:	COOLING S	YSTEM			Fluid Manu	Eluid HE		E ANTIER	
Unit Serial: UnivC Unit Worksite: Houston , TX Component Serial: UnivC Component Type: COOLING SYSTEM Fluid Grade: Interance for Lab No:: 190729012 Utated By: ENELSON ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dos of SCA. Resample at normal interval. ECTROCHEMICAL ANALYSIS IN PARTS PER MILLION Elements Ass SAMPLE SS Option 1 of SCA. Resample at normal interval. Interval. Status of SCA. Resample at normal interval.	Unit Seria: UnivC Unit Worksite: Houston, TX Fluid Grade: Fluid Grade: Enterson Fluid Grade: Fluid Grade: Component Seria: UniXC Component Seria: UniXC Component Seria: UniXC Fluid Grade: Fluid Grade: Component Seria: UniXC Sample Sample Sample One Seria: UniX Sample		Unit Model	: 3516				Component	onent Model:	3516				Brand/	Product:	WI DOTTIN		
Unit Worksit: Houston , TX Component Type: COOLING SYSTEM Intenance for Lab No.: 1907290612 ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT POAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dos of SCA. Resample at normal interval. CCROCHEMICAL ANALYSIS IN PARTS PER MILLION Eloments Res SAMPLE of of 00 DRAWN of of of 122 07/10/2019 c1 c1 c1 c1 131 06/28/2019 c1 c1 c1 c1 141 05/17/2019 c1 c1 c1 c1 142 02/27/2019 c1 c1 c1 c1 142 02/27/2019 c1 c1 c1 c1 143 05/17/2019 c1 c1 c1 c1 120 04/11/2019 c1 c1 c1 c1 121 07/10/2019 c1 c1 c1 c1 120 04/11/2019 c1 c1 c1 c1 121 05/17/2019 c1 c1 c1 c1 120 04/11/2019 <td>Unit Worksite: Houston, TX Component Type: COOLING SYSTEM internance for Lab No:: 10907290612 ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED! Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose of SCA. Resample at normal interval. CETROCHEMICAL ANALYSIS IN PARTS PER MILLION AB SAMPLE DRAWN db db</td> <td></td> <td>Unit Serial</td> <td>: UnivC</td> <td></td> <td></td> <td></td> <td>Comp</td> <td>onent Serial:</td> <td>UnivC</td> <td>122227</td> <td></td> <td></td> <td>Fluid</td> <td>Grade:</td> <td></td> <td></td>	Unit Worksite: Houston, TX Component Type: COOLING SYSTEM internance for Lab No:: 10907290612 ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED! Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose of SCA. Resample at normal interval. CETROCHEMICAL ANALYSIS IN PARTS PER MILLION AB SAMPLE DRAWN db		Unit Serial	: UnivC				Comp	onent Serial:	UnivC	122227			Fluid	Grade:			
ARALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dos of SCA. Resample at normal interval. CTROCHEMICAL ANALYSIS IN PARTS PER MILLION CTROCHEMICAL ANALYSIS IN PARTS PER MILLION	ANALYSIS INDICATES COOLANT MAINTENANCE/ SERVICE IS REQUIRED Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT POAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose of SCA. Resample at normal interval. ECTROCHEMICAL ANALYSIS IN PARTS PER MILLION Elements AB SAMPLE of some protection levels are LOW. Note SLIGHT DAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose of SCA. Resample at normal interval. CETROCHEMICAL ANALYSIS IN PARTS PER MILLION Elements AB SAMPLE of some protection levels are LOW. Note SLIGHT AB SAMPLE SAMPLE of some protection levels are LOW. Note SLIGHT AB SAMPLE of some protection levels are LOW. Note SLIGHT SAMPLE of some protection levels are LOW. Note SLIGHT SAMPLE of some protection levels are LOW. Note SLIGHT SAMPLE of some protection levels are LOW. Note SLIGHT SAMPLE of some protectine loss of heath anote some protection levels are LOW. Note SLIGHT																	
AB SAMPLE AB Joseph Joseph <thjoseph< th=""> Joseph Joseph</thjoseph<>	AB SAMPLE AP AP <th< th=""><th>SPECTR</th><th></th><th></th><th>RTS PER M</th><th>le at norm</th><th>al interval.</th><th></th><th></th><th>s, Note a cha</th><th>nge in the co</th><th>olant color. 1</th><th>i ne coolant a</th><th>ppears stable.</th><th>Suggest add</th><th>ding tuli main</th><th>tenance dose</th></th<>	SPECTR			RTS PER M	le at norm	al interval.			s, Note a cha	nge in the co	olant color. 1	i ne coolant a	ppears stable.	Suggest add	ding tuli main	tenance dose	
312 07/10/2019 <1 <1 <1 <1 <1 16 334 1619 58 207 65 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 </th <th>B12 07/10/2019 <1</th> <1	B12 07/10/2019 <1	SPECTR	OCHEMICAL ANA	LYSIS IN PA	ARTS PER M	le at norm	al interval.			Ele	ments	olant color. 1	i ne coolant a	ppears stable.	Suggest add	ding tuli main	tenance dose	
081 06/28/2019 <1	081 06/28/2019 <1	LAB NO.	OCHEMICAL ANA SAMPLE DRAWN		ARTS PER M	LLION	al interval.	Jan	Silloon	Ele	ments	unisseno	une coolant a	ppears stable.	Suggest add	ding tuli main	units and a second	
741 05/17/2019 <1	741 05/17/2019 <1	LAB NO. 0612	SAMPLE DRAWN 07/10/2019	LYSIS IN PA	ARTS PER M	LLION	al interval.	ی جا	400mg 16	Ele 6 334	ments	unissing SS	4000 1000 1000 1000 1000 1000 1000 1000	es	Suggest add	ung tuli main	^(u) see by <1	
320 04/11/2019 <1	820 04/11/2019 <1	LAB NO. 0612 1081	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019	LYSIS IN PA	ARTS PER M	LLION	al interval.	₹ <1	400 16 15	Ele 20 334 335	ments 	58 58	100 coolant a	es 65 68	Suggest add	ung tull main ung tull main solo tu tu tu tu tu tu tu tu tu tu tu tu tu	447773 447773 4467 447773 4467 447 447 447 447 447 447 447 447 44	
3/3 03/25/2019 <1 <1 <1 10 313 1330 42 204 01 <1 <1 <1 <1 <1 <1 16 313 1330 42 204 01 <1 <1 <1 <1 <1 <1 16 313 1330 42 204 01 <1 <1 <1 <1 <1 <1 16 341 1676 52 218 64 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	Solution Stand	LAB NO. 0612 1081 0741	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019	LYSIS IN P4 4 4 41 41 41 41 41	ARTS PER M	Cline at norm	al interval.	ي جا	400 116 115 117	Ele 6 334 335 359 259	ments 5 1619 1747	58 58 55 58	207 187 226	65 68 67	Suggest add	ung tuli main Ung tuli main Ung Ung Ung Ung Ung Ung Ung Ung Ung Un	Lenance dose	
MPLE INFORMATION Age UoM Filter Lub ab Sample Unit Lube UOM Filter Lub to: Drawn Time Age UOM Filter Lub til: 07/10/2019 34829 34829 HR No Ser til: 06/28/2019 34511 HR No Ser til: 05/17/2019 33541 HR No Ser til: 06/28/2019 34511 HR No Ser til: 06/28/2019 33541 HR No Ser til: 06/28/2019 33541 HR No Ser til: 06/28/2019 33541 HR No Ser til: 04/11/2019 32674 HR No Ser til: 04/11/2019 32674 HR No Ser til: 04/11/2019 32674 HR No Ser	MPLE INFORMATION Lub Sample Unit Lube UOM Filter Chgd. Ser No. Drawn Time Age UM Filter Chgd. Ser Ser 102 07/10/2019 34829 34829 HR No 5 103/28/2019 34541 HR No 5 103/28/2019 32574 HR No 5 103/28/2019 32272 HR No 5 103/28/2019 32272 HR No 5 103/28/2019 3277 HR NO 5 10	LAB NO. 0612 1081 0741 0820 0922	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 02/26/2010	(1) (1) (1) (1) (1) (1)	ARTS PER M	LLION	al interval.	ي جا	40 16 15 17 17	Ele 234 334 335 359 360 212	ments 1619 1659 1747 1752	58 58 55 59	207 187 228 208	65 68 67 66	Suggest add		 41 4	
Ab Sample Drawn Unit Time Lube Age UOM Filter Chgd. Lub Ser Ser Ser <td>Lab Sample Unit Lube UOM Filter Cup No. Sample Drawn Time Age UOM Filter Cup No. Sample Drawn Time Age UOM Filter Cup Sample Origonal 06/28/2019 34829 HR No Ser <t< td=""><td>LAB NO. 0612 1081 0741 0820 0873 0542</td><td>OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 03/25/2019 02/27/2019</td><td>LYSIS IN PA</td><td>ARTS PER M CA. Resamp ARTS PER M CI CI CI CI CI CI CI CI CI CI</td><td>LLION</td><td>al interval.</td><td><i>₹</i> <1</td><td>400 16 15 17 17 16 16</td><td>Ele 334 335 359 360 313 341</td><td>ments 1619 1659 1747 1752 1536 1676</td><td>58 58 58 59 42 52</td><td>207 187 226 228 204 218</td><td>65 65 66 67 66 61 64</td><td>Suggest add</td><td></td><td>41 41 41 41 41 41 41 41 41 41</td></t<></td>	Lab Sample Unit Lube UOM Filter Cup No. Sample Drawn Time Age UOM Filter Cup No. Sample Drawn Time Age UOM Filter Cup Sample Origonal 06/28/2019 34829 HR No Ser Ser <t< td=""><td>LAB NO. 0612 1081 0741 0820 0873 0542</td><td>OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 03/25/2019 02/27/2019</td><td>LYSIS IN PA</td><td>ARTS PER M CA. Resamp ARTS PER M CI CI CI CI CI CI CI CI CI CI</td><td>LLION</td><td>al interval.</td><td><i>₹</i> <1</td><td>400 16 15 17 17 16 16</td><td>Ele 334 335 359 360 313 341</td><td>ments 1619 1659 1747 1752 1536 1676</td><td>58 58 58 59 42 52</td><td>207 187 226 228 204 218</td><td>65 65 66 67 66 61 64</td><td>Suggest add</td><td></td><td>41 41 41 41 41 41 41 41 41 41</td></t<>	LAB NO. 0612 1081 0741 0820 0873 0542	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 03/25/2019 02/27/2019	LYSIS IN PA	ARTS PER M CA. Resamp ARTS PER M CI CI CI CI CI CI CI CI CI CI	LLION	al interval.	<i>₹</i> <1	400 16 15 17 17 16 16	Ele 334 335 359 360 313 341	ments 1619 1659 1747 1752 1536 1676	58 58 58 59 42 52	207 187 226 228 204 218	65 65 66 67 66 61 64	Suggest add		41 41 41 41 41 41 41 41 41 41	
No. Drawn Time Age Chgd. ser Sil2 07/10/2019 34829 34829 HR No S No1 06/28/2019 34511 HR No S V41 05/17/2019 33541 HR No S 120 04/11/2019 32674 HR No S 120 04/11/2019 32674 HR No S 120 04/11/2019 32674 HR No S	No.DrawnTimeAgeChgd.Ser61207/10/20193482934829HRNoSe08106/28/201934511HRNoSe74105/17/201933541HRNoSe82004/11/201932674HRNoSe87303/25/201932272HRNoSe67403/25/201932272HRNoSe87303/25/201931646HRNoSe	LAB NO. 0612 1081 0741 0820 0873 0542	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 INEORMATION	LYSIS IN PA	ARTS PER M UNIT VIIII VIIIII VIIII VIIII VIIII VIIIII VIIIII VIIII VIIII VIIII VIIII VIIIII VIIII VIIIII VIIIII VIIIII VIIIII VIIIIII VIIIII VIIIIIIII	e at norm	al interval.	<i>₹</i> <1	40 16 15 17 17 16 16	Ele 5 5 5 5 5 5 5 5 5 5 5 5 5	ments 1619 1659 1747 1752 1536 1676	58 58 58 59 42 52	207 226 228 204 218	65 68 67 66 61 64	Suggest and	400 Tull main 400 Tull main 400 Tull main 41 41 41 41 41 41 41 41 41 41	41 41 41 41 41 41 41 41 41 41 41	
$\begin{array}{ccccccc} & & & & & & & & & & & & & & & &$	$\begin{array}{cccccccccccccccccccccccccccccccccccc$	LAB NO. 0612 1081 0741 0820 0873 0542 SAMPLE Lab	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 INFORMATION Sample	LYSIS IN PA	ARTS PER M UNITED VIEW CI CI CI CI CI CI CI CI CI CI	LLION	al interval.	ية <1	60 16 15 17 17 16 16 16	Ele	ments 1619 1659 1747 1752 1536 1676	58 58 58 65 59 42 52	207 187 226 228 204 218	65 68 67 66 61 64	Suggest add	4000 Tull main 4000 Tull 4000 Tull 400	4038 40798 40998 40998 40 41 41 41 41 41 41 41 41 41 41 41 41 41	
141 05/17/2019 33541 HR No s informe. Para imprimir, seleccione el icono de la	05/17/2019 33541 HR No informe. Para imprimir, seleccione el icono de la 06/11/2019 32674 HR No impresora o, utilizando el teclado, haga clic en el 07/27/2019 31546 HR No impresora o, utilizando el teclado, haga clic en el	LAB NO. 0612 1081 0741 0820 0873 0542 CAMPLE Lab NO. 0612	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 INFORMATION Sample Drawn 07(10/2019	LYSIS IN PA	ARTS PER M CA. Resamp ARTS PER M CI CI CI CI CI CI CI CI CI CI	LLION	al interval.	₹ <1	00 16 15 17 17 16 16 16 16	Ele 5 334 335 359 360 313 341 TIES 2 UIDA	ments 1619 1659 1747 1752 1536 1676	58 58 55 59 42 52	207 187 226 228 204 218	65 68 67 66 61 64	Suggest add <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	ang tul main ⁴⁰ 0000 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	Clipton Clipto	
120 04/11/2019 32674 HR No since a set of the standard of the	^{B20} 04/11/2019 32674 HR № S B73 03/25/2019 32272 HR № S 542 02/27/2019 31646 HB № S	LAB NO. 0612 1081 0741 0820 0873 0542 SAMPLE Lab No. 0612 1081	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 INFORMATION Sample Drawn 07/10/2019 06/28/2019	LYSIS IN PA	ARTS PER M 41 41 41 41 41 41 41 41 41 41	LLION LLION V V V V V U V M HR HR	al interval.	,₹ <1 Fu	60 16 15 17 17 16 16 16 16 16 16 20 20 20 20 20 20 20 20 20 20 20 20 20	Ele 5 5 5 5 5 5 5 5 5 5 5 5 5	ments 1619 1659 1747 1752 1536 1676 NUEV	58 58 55 59 42 52	207 197 226 228 204 218	65 68 67 66 61 64 CON 6	Suggest add <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	<pre>dung tull main</pre>	unperformance unperformance <1	
	ara 03/25/2019 32272 HR No s impresora o, utilizando el teclado, haga clic en el	LAB NO. 0612 1081 0741 0820 0873 0542 CAMPLE Lab No. 0612 1081 0741	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 05/17/2019 04/11/2019 03/25/2019 02/27/2019 INFORMATION Sample Drawn 07/10/2019 06/28/2019 05/27/2019	LYSIS IN PA	ARTS PER M 	LLION LLION V V V V UOM HR HR HR HR	al interval.	₹ <1 Se	00 16 15 17 17 16 16 16 16 16 17 17 16 16 16 17 17 17 16 16 16 17 17 17 16 16 15 17 17 17 17 16 16 15 17 17 17 16 16 15 17 17 17 16 16 15 17 17 16 16 15 17 17 17 16 16 16 15 17 17 16 16 16 16 17 17 16 16 16 16 17 17 16 16 16 16 16 16 17 17 16 16 16 16 16 16 16 16 17 17 16 16 16 16 16 16 16 16 16 16	Ele 5 334 335 359 360 313 341 TIES Á UNA Para	ments 1619 1659 1747 1752 1536 1676 NUEV	58 58 55 59 42 52	207 187 226 228 204 218 Staña	65 68 67 66 61 64 CON 6	Suggest and <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1 <1	ding full main	Cl C	
	1111/16501a 0, utilizando enteciado, naga circ en el	LAB NO. 0612 1081 0741 0820 0873 0542 AMPLE Lab No. 0612 1081 0741 0820	OCHEMICAL ANA SAMPLE DRAWN 07/10/2019 06/28/2019 04/11/2019 03/25/2019 02/27/2019 02/27/2019 02/27/2019 02/27/2019 06/28/2019 06/28/2019 06/28/2019 06/27/2019 04/11/2019	LYSIS IN PA	ARTS PER M 41 41 41 41 41 41 41 41 41 41	LLION LLION V V V V UOM HR HR HR HR HR	al interval.	,≰ <1 Se info	do do do do do rme.	Ele 334 335 359 360 313 341 TIES á una Para	ments 1619 1659 1747 1752 1536 1676 NUEV	58 58 59 42 52 70 70 70 70 70 70 70 70 70 70 70 70 70	207 197 226 228 204 218 staña selecci	65 68 67 66 61 64 CON 6 Cione	<pre></pre>	 cling full main cling cling cling <licling< li=""> cling cling clin</licling<>	Internance dose	

Para descargar, seleccione el icono de descarga.

This analysis is intended as an aid in predicting mechanical wear. Test results, ma and apply only to this sample as provided. No guarantee, expressed or implied, is equipment and all of its components is the responsibility of the equipment owner. Testing performed by Bureau Veritas, an ISO/IEC 17025-2017 accredited laboratory by A accreditation. Click here for Tests and Methodologies





RECOMMENDATIONS

Maintenance for Lab No.: 201907290612 Evaluated By: ENELSON ANALYSIS INDICATES COOLANT MAINTENANCE / SERVICE IS REQUIRED! Note some phosphorous present. Corrosion protection levels are LOW. Note SLIGHT FOAMING, which can cause loss of heat transfer if this level increases, Note a change in the coolant color. The coolant appears stable. Suggest adding full maintenance dose of SCA. Resample at normal interval.

> Para enviar un informe por correo electrónico desde la pantalla Detalles, pulse el botón *ENVIAR CORREO ELECTRÓNICO*.

ENVIAR UN INFORME POR CORREO ELECTRÓNICO



Zinc

horus

CANCEL

SEND

Calcium

Magnesiun

	IECK		-	Ip ⑦ Search Search
Transform Transform	Saca may gance		EMAIL SAMPLE REPORT	×
HOME ACT	IONS SAMPLES LABELS	EQUIPMENT AN	Recipient(s) Email*	
← ВАСК	K TREND GRAPH	SEND EMAIL	Type recipient email and press Enter or search contacts	Sample Condition Status 🧲
			Subject*	
	Lab Custaman ID#	210102	Oil Analysis Report - ACME Demo - Houston , TX - 1107/ELMER FUDD EAST/MAIN - COOLING SYSTEM	
	Lab Customer ID#: 319182 Company Name: ACME Demo		Message*	Nork Order No.:
	Address:	123,		Filter Age: 0
		Editiontion, Alber	I,	22
	UNIT INF	ORMATION	Acme Demo has e-mailed you a copy of an oil analysis report. Please see the attached pdf report.	FLUID INFORMATION
	Unit ID: Unit Manufacturer:	1107/ELMER FUE Caterpillar		d Manufacturer: CHEVRON Brand/Product: HEAVY DUTY PF ANTIFR
	Unit Model: Unit Serial:	3516 UnivC		Fluid Grade:
	Unit Worksite:	Houston , TX		
PECONNE	NEATIONS			
Maintenan	ice for Lab No.:	ANALYSIS INDICA		. Note SLIGHT FOAMING, which can cause loss of heat
201907290 Evaluated	0612 By: ENELSON	transfer if this leve		ample at normal interval.
Results	History Actions	Graphs Atta	To reply please contact: acme_demo@bureauveritas.com	
COLOR			Thank you,	
SPECI	So obrirá u		a ventana. Haz elie en el icena	
	Se abilia u		a ventana. Haz ciic en el icono	

Buscar para encontrar y añadir destinatarios desde tu Agenda o escribe el correo electrónico del destinatario y haz clic en Intro. Añade un mensaje y pulsa el botón **ENVIAR**.



SERVICIO DE SOPORTE DE CASTROL LABCHECK: Teléfono: 866-LABCHECK (522-2432) Labchecksupport@bureauveritas.com https://www.labcheckresources.com/



